Unpacking, Assembly & Installation

Unpacking

Inspect the delivered box for any external sign of damage. If any is found make a note of this before opening the box.

Remove all parts in the delivery and check each item against the delivery note. Note any parts that may be missing or deviations from the packing list.

Inspect all parts for any sign of damage and if any are found make a note of all damage.

Any missing parts or damage must be reported to the supplier and delivery agent as soon as possible.

Assembly

The top side of the AutoQuest autosampler, sample carousel platen and needle holder are shown in Figure 1

- a) Place the sample carousel platen onto the drive spindle ensuring that the locating pin is correctly positioned and located.
- b) Remove the holding nut and washer from the needle assembly and put the needle (caution: sharp needle) through the hole of the needle holder and replace the nut and washer and tighten the nut so that the needle is fixed firmly in place.
- c) The sample feed tube from the needle to the Rheodyne valve is terminated with a short length of FEP tube onto which is swaged a stainless ferrule. This is placed in port No. 5 of the Rheodyne valve and the captive nut is screwed into this port and tightened to give a tight, leak free joint.

Figure 1

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